


A B C D E F G H J

1  
2  
3  
4  
5

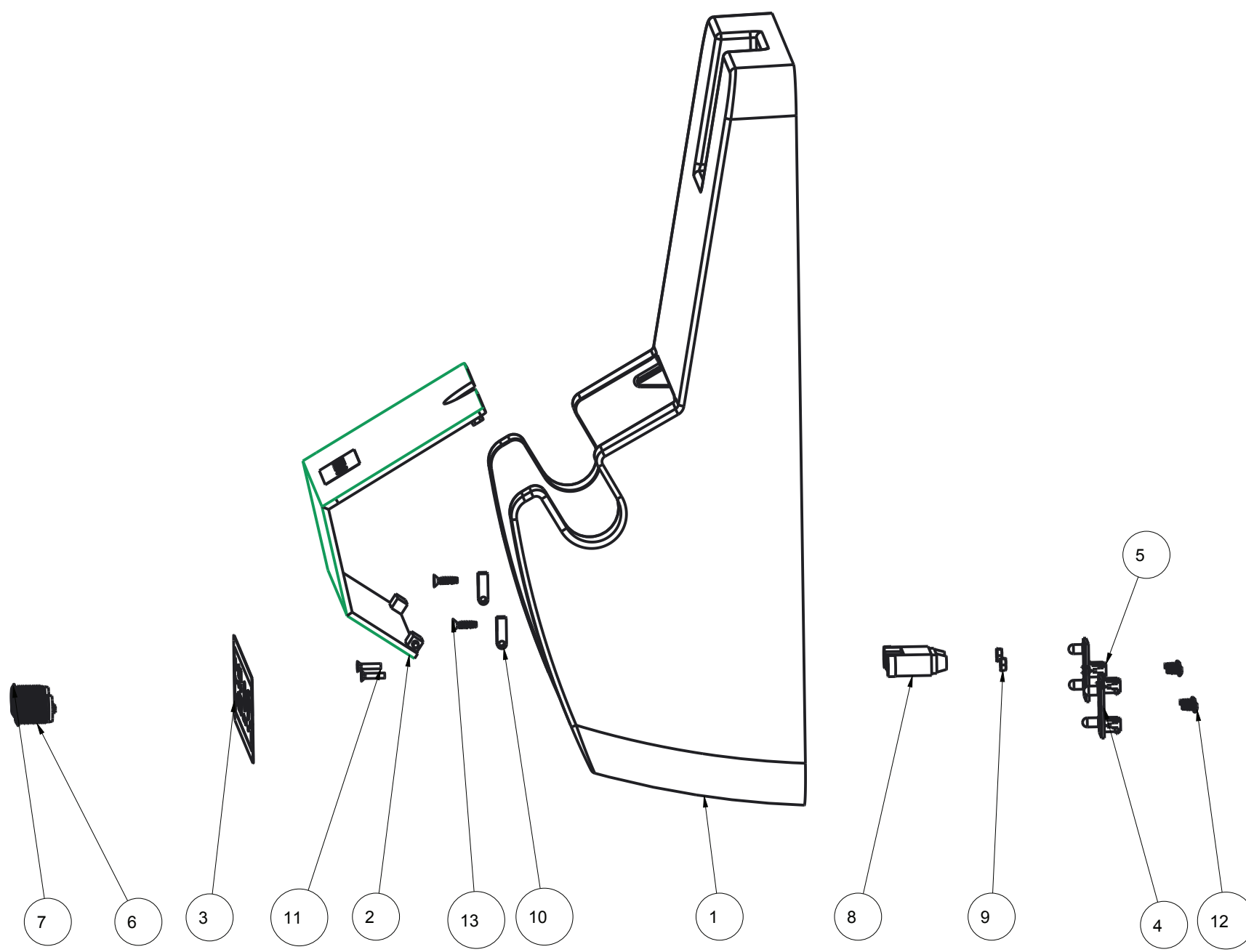

13	5510-0142	THREAD FORMING TORX FLAT HEAD 6-19 X 1/2	2
12	5510-0141	THREAD FORMING PAN HEAD 4-40 X 1/4	4
11	5510-0033	S/S PFH SCREW #4-40 X 1/2" LG	2
10	3200-0250	HINGE BLOCK, COVER	2
9	5535-0006	#4-40 KEPS NUT, S.S	2
8	2700-0037	FRONT PANEL USB CABLE	1
7	2700-0035	RELAY POWER SWITCH	1
6	2700-0036	COMPUTER POWER SWITCH	1
5	5900-0078	INDICATOR BOARD, TEMPERATURE	1
4	5900-0077	INDICATOR BOARD, POWER	1
3	7401-0009	OVERLAY, FRONT	1
2	3250-0136	COVER, FRONT	1
1	3350-0009	BEZEL, FRONT	1
ITEM	PART NO.	DESCRIPTION	QTY

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO OTHERS FOR ANY PURPOSE NOR USED FOR MANUFACTURING PURPOSES WITHOUT WRITTEN PERMISSION FROM DISTEK.

**DISTEK**  
121 NO. CENTER DR.  
NORTH BRUNSWICK, N.J. 08902

TITLE		SCALE	
FRONT BEZEL ASSEMBLY		1/1	
OPT-DISS 410		SIZE	SHEET OF REV
		B	11 A
		DRAWING NO.	
		2800-0491	