
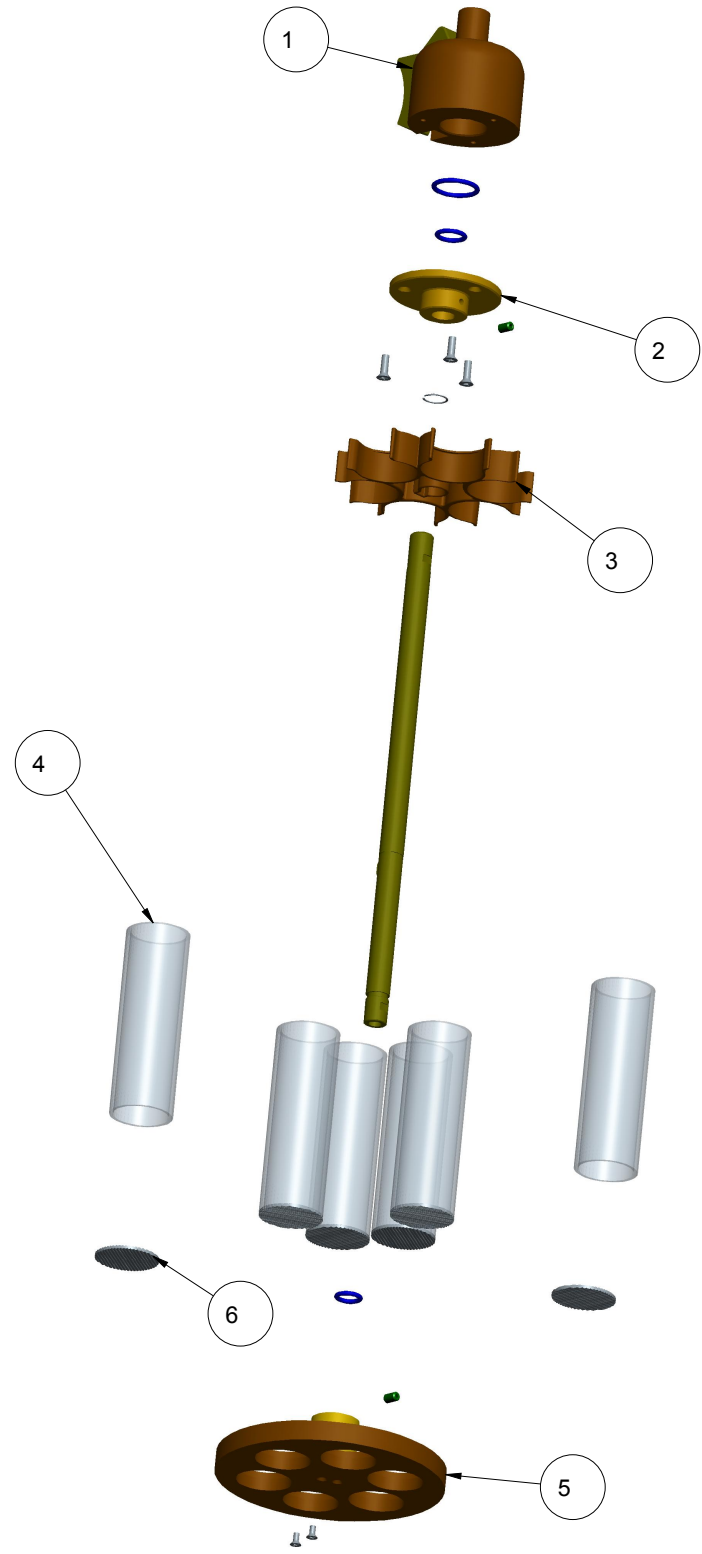


A B C D E F G H J

AUTO DETECT BASKET ASSY 2800-0459		
ITEM	PART NUMBER	REPLACE WITH
1	2800-0462	2800-0466
5	2800-0467	2800-0468

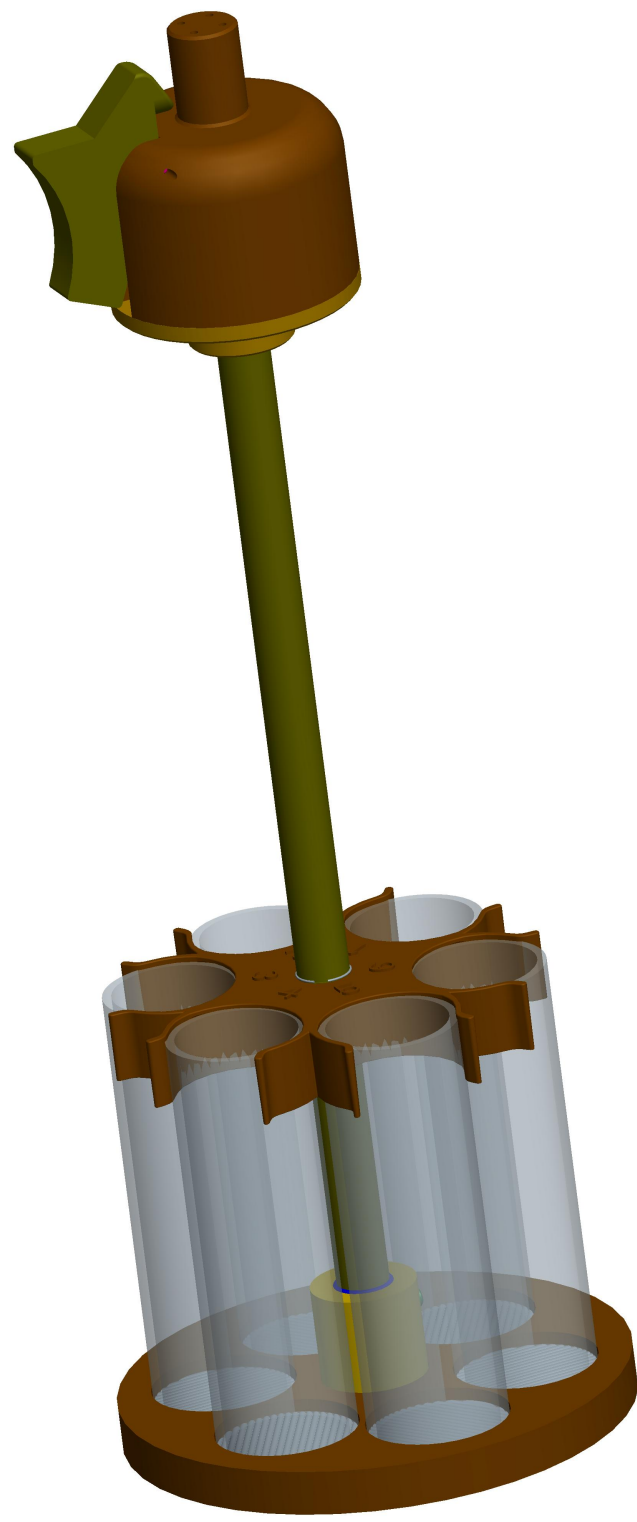

ITEM	PART NO.	DESCRIPTION	QTY
6	3200-0032	BASKET SCREEN 10 MESH	6
5	2800-0467	BOTTOM FLANGE ASSY, STD	1
4	3010-0024	BASKET TUBE	6
3	3350-0375	TOP FLANGE, BASKET	1
2	3200-0301	BASKET CONNECTOR COVER	1
1	2800-0462	BASKET CONNECTOR ASSY, STD	1

CONFIDENTIAL INFORMATION  
 THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION  
 AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO  
 OTHERS FOR ANY PURPOSE NOR USED FOR  
 MANUFACTURING PURPOSES WITHOUT WRITTEN  
 PERMISSION FROM DISTEK.

**DISTEK**  
 121 NO. CENTER DR.  
 NORTH BRUNSWICK, N.J. 08902

TITLE  
**BASKET ASSEMBLY**  
 sensIR 3200, STD

SCALE **1/2**  
 SIZE **B** SHEET OF **11** REV **D**  
**REVIEW ONLY**

1  
2  
3  
4  
5