
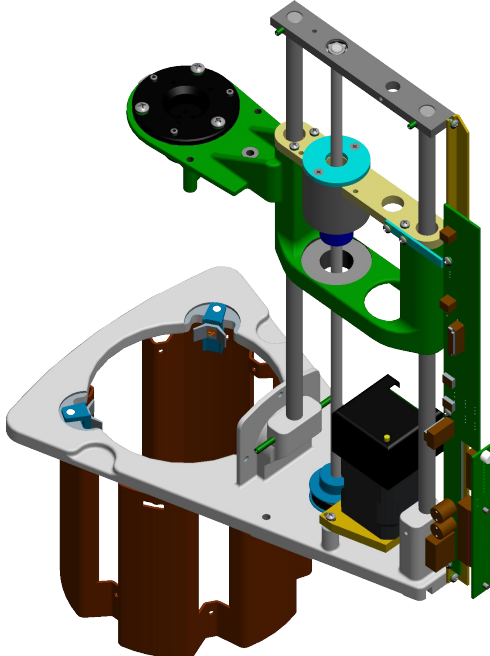


A B C D E F G H J

MODULE ASSEMBLIES FOR VOLTAGES AND DETECTION				
ITEM	PART NUMBER	115V, AUTO	230V, STD	230V, AUTO
		2800-0470	2800-0456	2800-0474
REPLACE WITH				
1	2800-0461	2800-0471	2800-0461	2800-0471
2	2600-0329	2600-0329	2600-0331	2600-0331
4	2800-0464	2800-0475	2800-0464	2800-0475
6	2800-0463	2800-0476	2800-0463	2700-0476

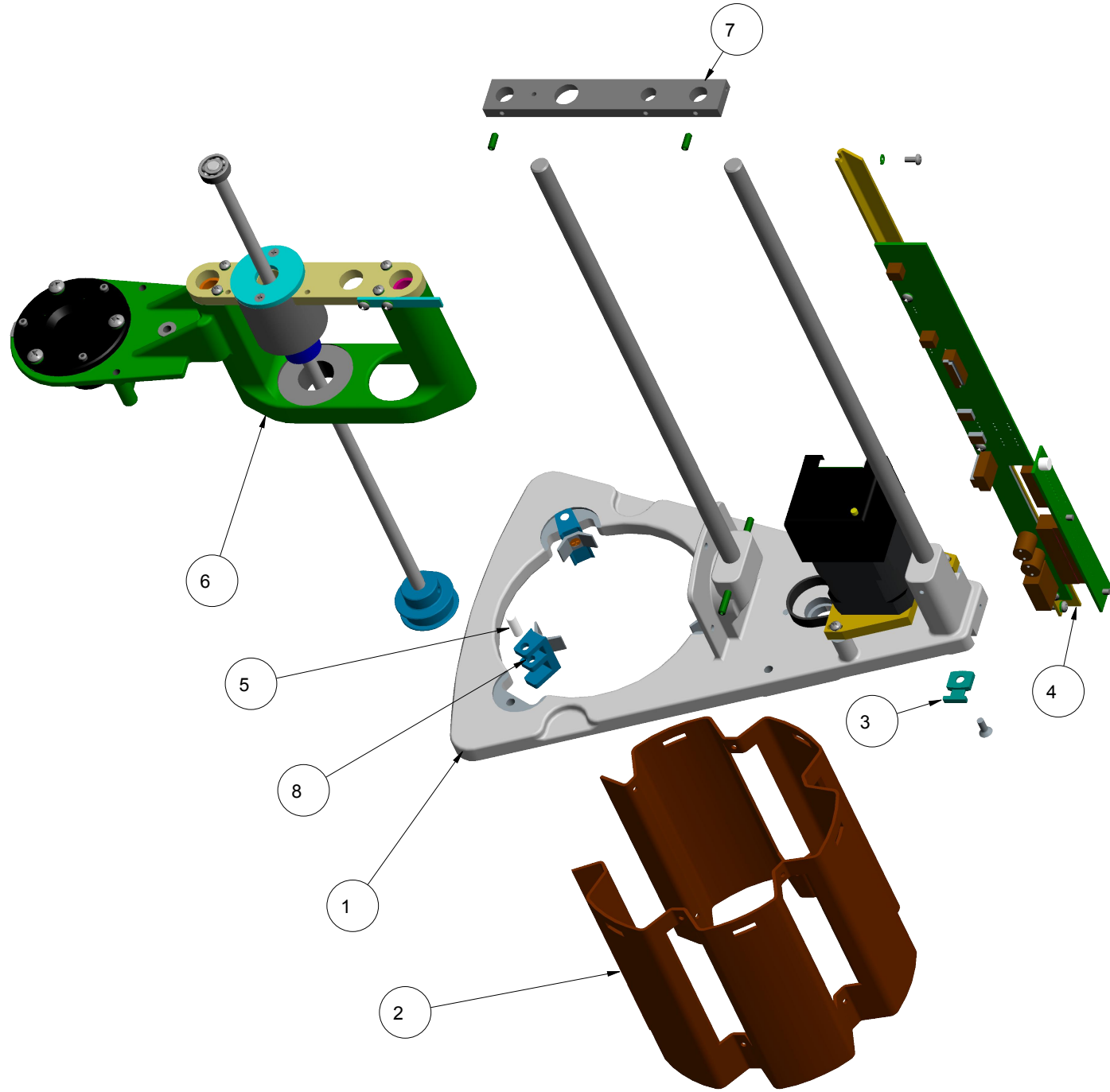

8	2851-0016	VESSEL CENTERING ARM ASSEMBLY	3
7	3200-0220	TOP PLATE	1
6	2800-0463	STROKE SYSTEM ASSEMBLY	1
5	5545-6068	PIN, CENTERING ARM, 3/4"	3
4	2800-0464	PCB ASSEMBLY	1
3	5570-1002	NYLON CABLE HOLDER	2
2	2600-0329	HEATER ENCLOSURE ASSEMBLY, 115V	1
1	2800-0461	DRIVE SYSTEM ASSEMBLY	1
ITEM	PART NO.	DESCRIPTION	QTY

<p>CONFIDENTIAL INFORMATION THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO OTHERS FOR ANY PURPOSE NOR USED FOR MANUFACTURING PURPOSES WITHOUT WRITTEN PERMISSION FROM DISTEK.</p>		<p>121 NO. CENTER DR. NORTH BRUNSWICK, N.J. 08902</p>	
TITLE		SCALE	
<p>MODULE ASSY, 115V, STD sensIR 3200</p>		<p>1/3</p>	
		SIZE	SHEET OF REV
		B	11 D
REVIEW ONLY			

1  
2  
3  
4  
5