

A B C D E F G H J


1

2

3


4

5

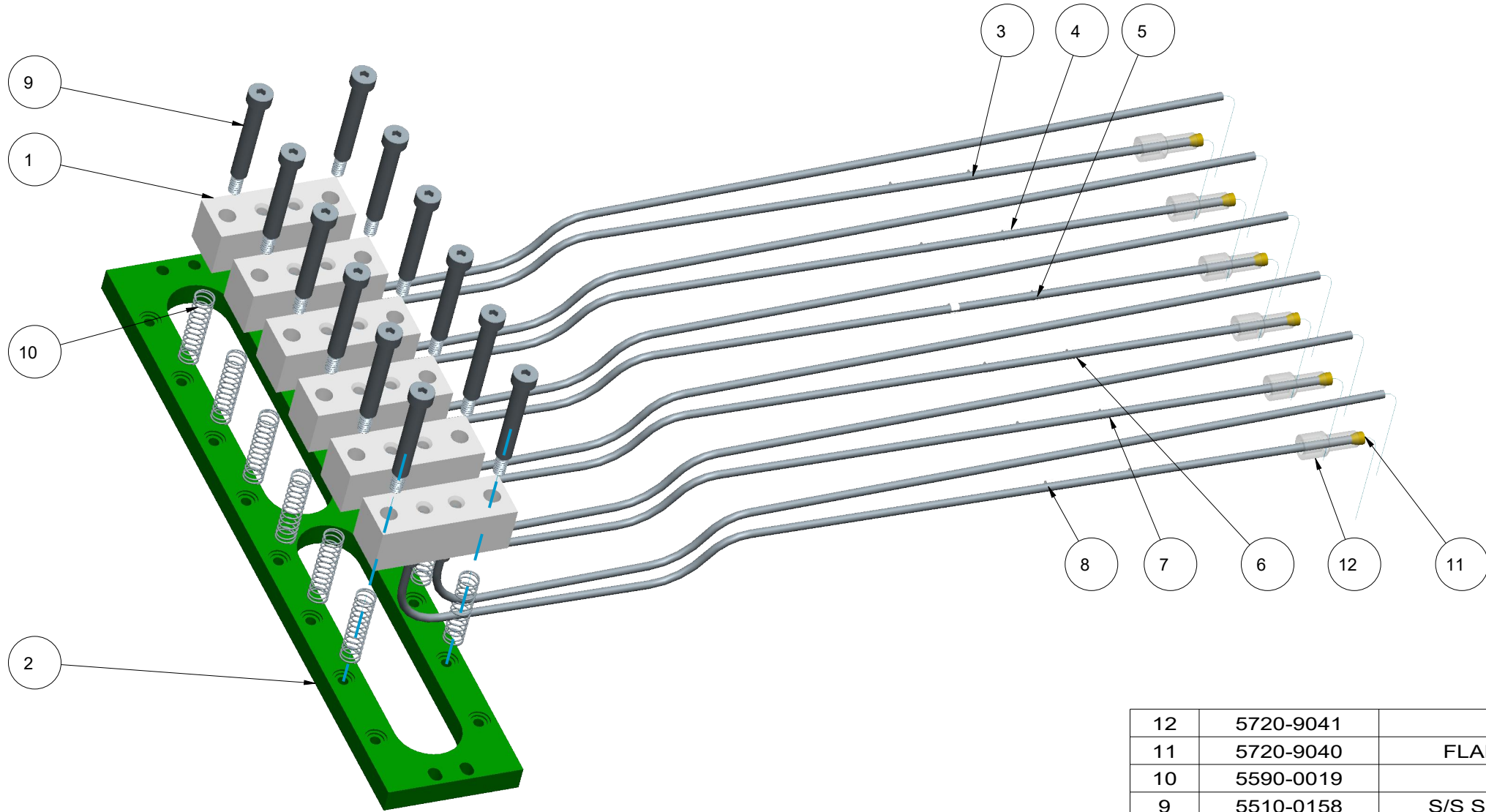

POSITION	SLEEVE COLOR
1	RED
2	WHITE
3	YELLOW
4	GREEN
5	BLUE
6	BLACK
7	CLEAR
8	NONE

12	5720-9041	1/4-28 FITTING, 1/8 OD TUBING	6
11	5720-9040	FLANGELESS FERRULE, 1/8 OD TUBING	6
10	5590-0019	SPRING, DIST. BLOCK MOUNT	12
9	5510-0158	S/S SHOULDER SCREW, 1/4 DIA. x 1-1/2 Lg.	12
8	4958-0003-6	TUBING MARKING, SNAP ON, #6	2
7	4958-0003-5	TUBING MARKING, SNAP ON, #5	2
6	4958-0003-4	TUBING MARKING, SNAP ON, #4	2
5	4958-0003-3	TUBING MARKING, SNAP ON, #3	2
4	4958-0003-2	TUBING MARKING, SNAP ON, #2	2
3	4958-0003-1	TUBING MARKING, SNAP ON, #1	2
2	3200-0313	DISTRIBUTION PLATE	1
1	2800-0383	DISTRIBUTION BLOCK ASS'Y	6
ITEM	PART NO.	DESCRIPTION	QTY

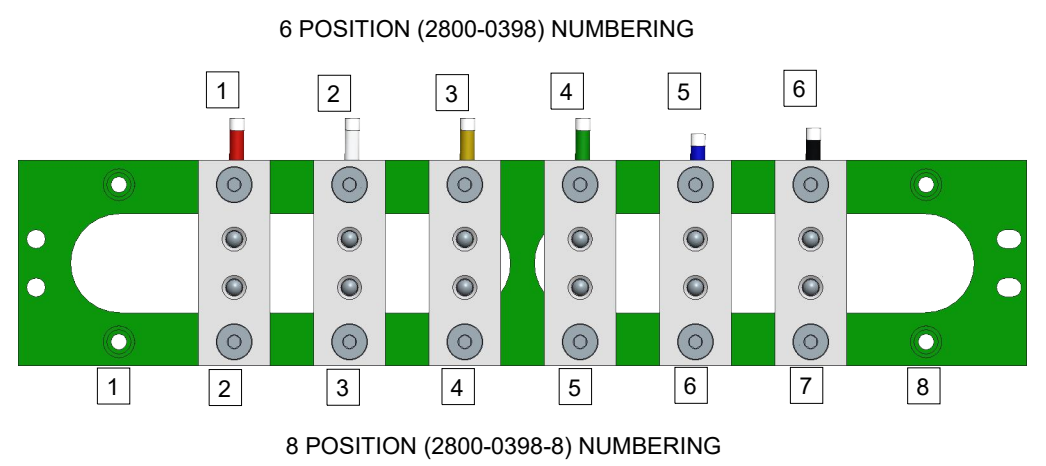

CONFIDENTIAL INFORMATION
 THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION AND SUCH INFORMATION IS NOT TO BE DISCLOSED TO OTHERS FOR ANY PURPOSE NOR USED FOR MANUFACTURING PURPOSES WITHOUT WRITTEN PERMISSION FROM DISTEK.

DISTEK
 121 NO. CENTER DR.
 NORTH BRUNSWICK, N.J. 08902

TITLE: **SAMPLE DISTRIBUTION ASS'Y**
ECLIPSE 5300

SCALE: **1/2**

SIZE: **B** SHEET OF **11** REV. **C**

REVIEW ONLY