

Date	7 Oct 2011	Service Note #	7100-1015	Updated	
Product	7100	Description	Removing Module		
Created By	Andrew Infante				

Release	<input checked="" type="checkbox"/>	Internal	<input checked="" type="checkbox"/>	Distributors	<input type="checkbox"/>	Customers
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	--------------------------	------------------

Parts Required	Serial Numbers Affected
Phillips Head Screw Driver	N/A

Steps

- 1 – Remove the vessel cover of the module you wish to remove.

- 2 – Raise shaft of the module.


- 3 – Power down unit

- 4 – Remove vessel.


- 5 –Using a Phillips head screw driver remove the three screws above the vessel plate.




5 – Remove top screw.

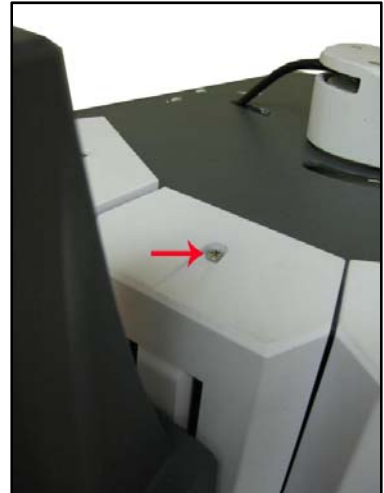

6 – Gently pull height adjustment lever and slide cover up.

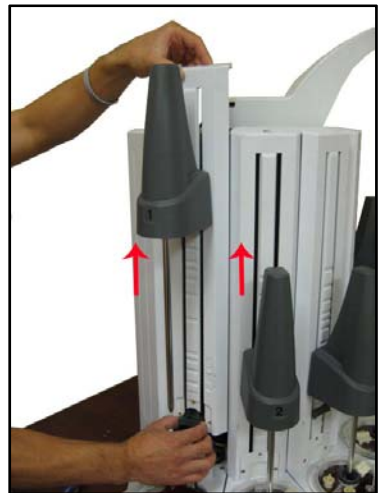

7 – Remove two screws from vessel plate.

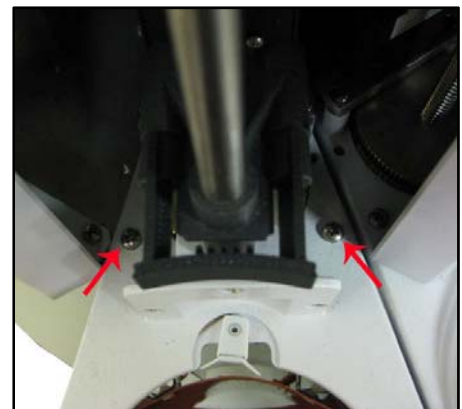

8 – Using vessel plate and guide rod gently pull the position straight out.

