

| | | | | | |
|-------------------|----------------|-----------------------|---------------------|----------------|--|
| Date | 7 Oct 2011 | Service Note # | 7100-1016 | Updated | |
| Product | 7100 | Description | USB Firmware Update | | |
| Created By | Andrew Infante | | | | |

| | | | | | | |
|----------------|-------------------------------------|-----------------|-------------------------------------|---------------------|--------------------------|------------------|
| Release | <input checked="" type="checkbox"/> | Internal | <input checked="" type="checkbox"/> | Distributors | <input type="checkbox"/> | Customers |
|----------------|-------------------------------------|-----------------|-------------------------------------|---------------------|--------------------------|------------------|

| | |
|-----------------------|--------------------------------|
| Parts Required | Serial Numbers Affected |
| USB Key | N/A |

Steps

1 – Firmware Upgrades will be sent from Distek via email. Once received, copy the file onto a USB Flashdrive.


2 – Insert USB Flashdrive into unit.



3 – Using the touch screen user interface navigate to Service by selecting:

Settings>System Settings>Service

4 – Select the appropriate icon to upgrade as specified in Distek’s Firmware Release.

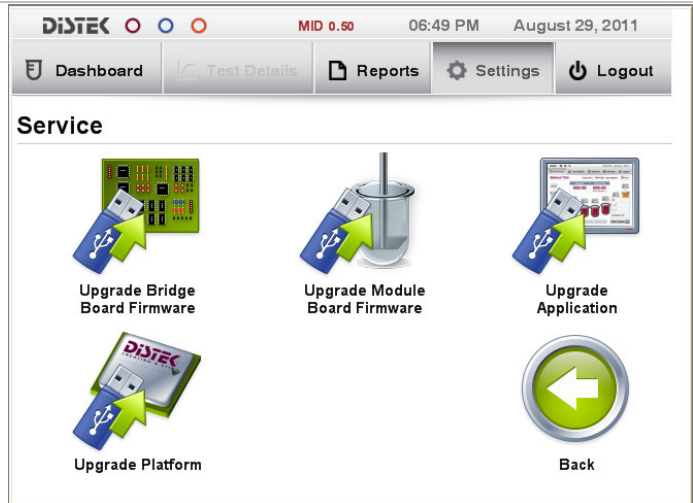

5 – When prompted enter password “symfw”.

6 – System will prompt that information has been found. Select continue. The files will automatically download and the system will reboot to complete the Firmware Upgrade.