


<b>Date</b>	05 Apr 2011		<b>Release Note #</b>	4300-1003	<b>ECO Ref #</b>	0847
<b>Product</b>	4300		<b>Description</b>	3-Way Distributor		
<b>Revision Change</b>	<b>From</b>	F	<b>To</b>	G	<b>Created By</b>	Joey Yangco

<b>Release</b>	<input checked="" type="checkbox"/>	<b>Internal</b>	<input checked="" type="checkbox"/>	<b>Distributors</b>	<input checked="" type="checkbox"/>	<b>Customers</b>
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

<b>Assembly / Part #</b>	2800-0127 Rev. G/ 3250-0041 Rev. G
<b>Serial #</b>	SP02504 or later
<b>Description</b>	Replace the 3-way syringe pump distributor from 3 pieces to one piece design.

This part was redesigned to incorporate all 3 existing parts (3-way diverter, adapter and aluminum piece) into a one piece design to simplify the assembly and increase reliability.

**Existing 3-Way Diverter**

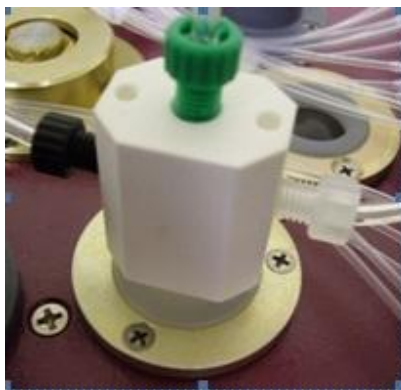

**One Piece 3-Way Diverter**





For older models using flat head screws to secure the 3-way distributor, three 4/40 x 0.375 inch long socket head screws will be required for replacement.

When using the one piece 3-way diverter on existing instruments, be aware to position the diverter in a way that you can access all the fittings without any inference from the adjoining diverters and the top cover.

---