

Date	01 Jul 2011		Release Note #	6x00-1007	ECO Ref #	0873
Product	Evolution 6x00		Description	pc/104 Firmware 2.17		
Revision Change	From	2.16	To	2.17	Created By	JSB

Release	<input checked="" type="checkbox"/>	Internal	<input checked="" type="checkbox"/>	Distributors	<input checked="" type="checkbox"/>	Customers
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

Assembly / Part #	9200-6163 (Firmware Revision)
Serial #	<p>Evolution 6100 (Bathless) – 6102056 or Later Evolution 6300 (Water Bath) – None</p> <p><i>Note: Due to a critical defect which was quickly detected only a few instruments utilized this release. Firmware 2.18 was released shortly after this version. This version is now obsolete and should no longer be utilized.</i></p>
Description	<p>In order to be compatible with the Distek Dosage Delivery System (DDS) the Firmware was updated with the following changes.</p> <ul style="list-style-type: none"> • The option “Tablet Dropper” was added to the “System Setup” screen. Which can be set to On or Off. • When set to On the user will be prompted during a Method run to introduce the dosage at the defined Temperature (once pre-heating is completed) or at a specified Time. <p>The Evolution 4300 and Cipher Software are compatible with this new Firmware and the Dosage Delivery System (DDS).</p>