



Release Note

Date	25 Jul 2011		Release Note #	6x00-1009	ECO Ref #	0885
Product	Evolution 6x00		Description	pc/104 Firmware 2.18		
Revision Change	From	2.17	To	2.18	Created By	JSB

Release	<input checked="" type="checkbox"/>	Internal	<input checked="" type="checkbox"/>	Distributors	<input checked="" type="checkbox"/>	Customers
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

Assembly / Part #	9200-6163 (Firmware Revision)
Serial #	Evolution 6100 (Bathless) – 6102083 or Later Evolution 6300 (Water Bath) – 6301399 or Later
Description	<ul style="list-style-type: none"> When using the Time mode function of the Evolution 6300 and the Dosage Delivery System, the Time screen was not displayed properly on the instrument. This defect is specific to the Evolution 6300 and is resolved in this Firmware release. In order to minimize the potential for the 6100 heaters to energize when the instrument is idle, the instrument sends a command every one second to ensure the heaters (triacs) are turned off. This is a preventative measure in rare cases where the DC 5.0V falls below the minimum threshold. This defect is specific to the Evolution 6100.