

Date	11 Jan 2012	Service Note #	MISC-1047	Revision	N/A
Product	2500x symphony 7100	Description	Networked Printing Limitations		
Created By	J. Yangco				

Release	<input checked="" type="checkbox"/>	Internal	<input checked="" type="checkbox"/>	Distributors	<input checked="" type="checkbox"/>	Customers
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

Parts Required	Serial Numbers Affected
N/A	2500x with Firmware Rev. 1.04 and older 7100 with Firmware Rev. 1.05 and older

Networked Printing Limitations

The following item describes the limitations when setting up the instrument to print using printer server:

1. IP settings are assigned automatically using DHCP only
2. Instrument must use the "Share Name" of the printer server
3. In the "Advanced Setup" screen, "Network Type" must only be set to DHCP. The "TCP-IP" setting is not supported at this time.