


Date	05 Feb 2013	Service Note #	Opt-Diss-1069	Revision	None
Product	Opt-Diss 405	Description	Wrong Revision in About Screen		
Created By	J. Yangco				

Release	<input checked="" type="checkbox"/>	Internal	<input checked="" type="checkbox"/>	Distributors	<input checked="" type="checkbox"/>	Customers
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

Parts Required	Serial Numbers Affected
N/A	Instruments with Software Revision 1.18.1 Build 3583

Wrong Revision in About Screen


Defect: The software version number that is displayed in the “About” screen does not match the software installed to run the Opt-Diss 405. The screen displays “OPT-DISS Version 1.18.0 Build 3583”.

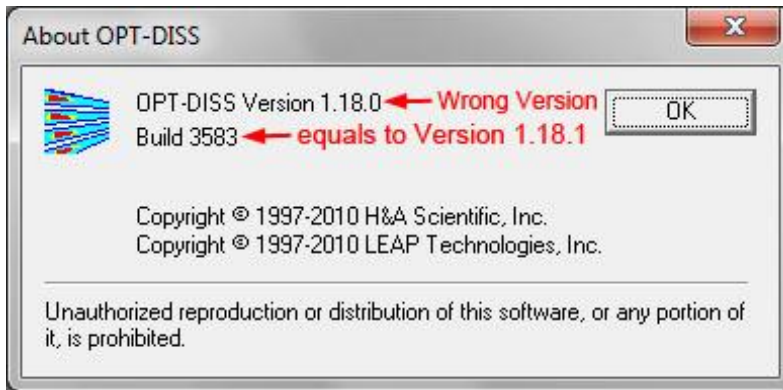

Opt-Diss “About” Screen Displays Wrong Software Version

Software Verification: The correct software revision can be verified using the “Build” number instead. The Opt-Diss 405 software Version 1.18.1 will have a build number of “3583”.

Another way to determine if the software is version is correct will be the Opt-Diss application executable date of 9/19/2011. You may be able to use these dates to determine the actual version of the executable, but these dates can change under certain circumstances.

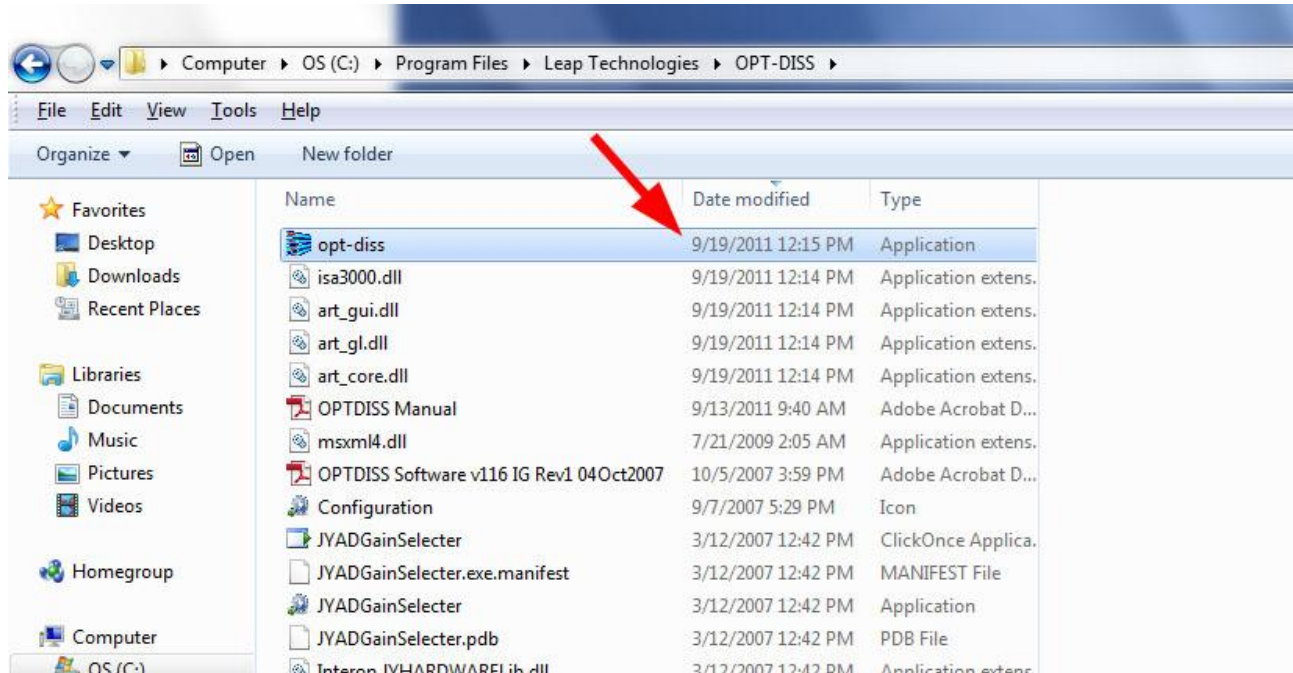

Opt-Diss Application Executable Date