

Date	04 Feb 2014	Service Note #	MISC-1095	Revision	NA
Product	Universal Basket Access Cap	Description	Universal Basket Access Cap O-Ring Installation		
Created By	J. Yangco				

Release	<input checked="" type="checkbox"/>	Internal	<input type="checkbox"/>	Distributors	<input type="checkbox"/>	Customers
----------------	-------------------------------------	-----------------	--------------------------	---------------------	--------------------------	------------------


Parts Required	Serial Numbers Affected
3250-0126 Rev. F 5570-8025 O-Ring 5570-8027 O-Ring	NA

Universal Cap O-Ring Installation

- Use current Universal Basket Cap, 3250-0126 Rev. F
- **2100, 6100 and 6300** instruments: use the side with the orange Silicone O-ring, 5570-8025 facing up
- **2500 and 7100** instruments: use the side with the thicker, black EPDM O-ring, 5570-8027 facing up
- Assembly with both O-rings installed: 2800-0055, Rev. D



2100, 6100 and 6300 Basket Access Cap Side

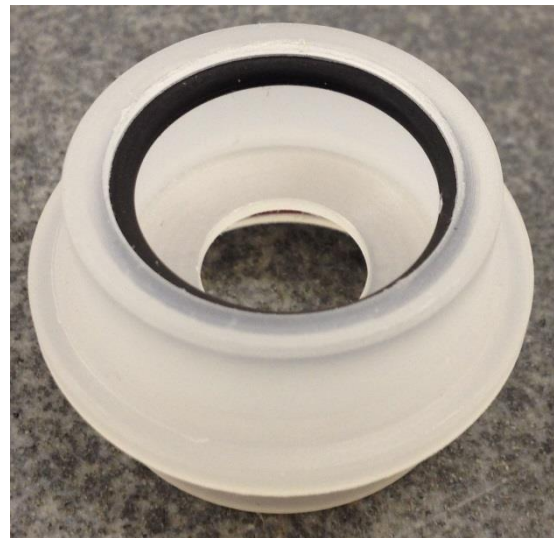

2500 and 7100 Basket Access Cap Side