

<b>Date</b>	05 May 2014	<b>Service Note #</b>	7100-1056	<b>Updated</b>	NA
<b>Product</b>	symphony 7100		<b>Created By</b>	Joey Yangco	
<b>Description</b>	Power Down Instrument When Removing and Installing Modules				

<b>Release</b>	<input checked="" type="checkbox"/>	<b>Internal</b>	<input checked="" type="checkbox"/>	<b>Distributors</b>	<input checked="" type="checkbox"/>	<b>Customers</b>
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

Parts Required	Serial Numbers Affected
None	All

## Power down instrument when removing or installing the module(s)

**Reason:** To avoid electrical surge on the module that can cause the module or instrument to malfunction.

**Pre-requisite:**

Service Engineer or person performing this service must have proper training in servicing the instrument.

**Solution / Action**



Before removing or installing the module(s), discharge any static electricity from your body by touching a bare screw from the instrument then turn off the instrument (you may also disconnect the AC power cord).