

| | | | | | |
|--------------------|--------------------------------|-----------------------|-------------------|----------------|----|
| Date | 18 May 2015 | Service Note # | Opt-Diss - 2045 | Updated | NA |
| Product | Opt-Diss 405 | | Created By | J. Yangco | |
| Description | Phase Time not getting updated | | | | |

| | | | | | | |
|----------------|-------------------------------------|-----------------|-------------------------------------|---------------------|-------------------------------------|------------------|
| Release | <input checked="" type="checkbox"/> | Internal | <input checked="" type="checkbox"/> | Distributors | <input checked="" type="checkbox"/> | Customers |
|----------------|-------------------------------------|-----------------|-------------------------------------|---------------------|-------------------------------------|------------------|

| Parts Required | Serial Numbers Affected |
|----------------|--|
| None. | Opt-Diss instrument with software Rev. 2.0.0 |

Phase time not getting updated

Reason:

The *Phase Time* in *Edit Sampling Time* dialog window is not updating, if the up/down buttons to change the # *Readings* or *Interval* are not used.

Solution / Action:

After manually entering the desired number of readings and interval time, the user must use the up/down buttons to have the software calculate and update the total phase time.