

<b>Date</b>	18 Jun 2015	<b>Service Note #</b>	OPT-DISS-2056	<b>Updated</b>	NA
<b>Product</b>	Opt-Diss 405		<b>Created By</b>	J. Yangco	
<b>Description</b>	3880-OPTDISS-OQ-S OPT-DISS Software OQ/PQ Validation Document (Rev. 3.9)				

<b>Release</b>	<input checked="" type="checkbox"/>	<b>Internal</b>	<input checked="" type="checkbox"/>	<b>Distributors</b>	<input checked="" type="checkbox"/>	<b>Customers</b>
----------------	-------------------------------------	-----------------	-------------------------------------	---------------------	-------------------------------------	------------------

<b>Parts Required</b>	<b>Serial Numbers Affected</b>
None.	All instruments that use software Rev. 2.0.0 and later


## Distek Dissolution Bath Configuration

### Reason:

When performing software validation, section 1.18 (File Menu), the requirements for File-Save As Method procedure does not report the bath settings if there is no Distek dissolution bath configured.

### Solution / Action

1. Login to the Opt-Diss software as part of the optra-administrators group.
2. On the main menu, select *Tools | System Configuration*. Configure the *Bath Configuration* (Bath 1 and Bath 2 if necessary) as shown:

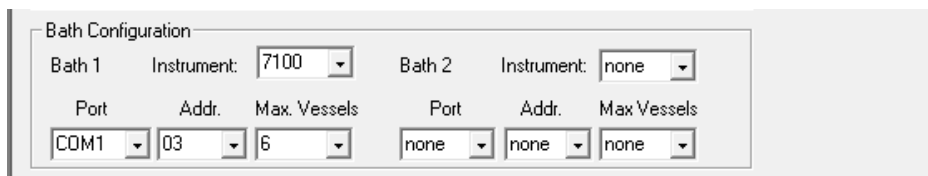

**Bath configuration setting**

3. Select OK and enter the password to accept the changes.



**After completing the validation, change the Bath Configuration to its original setting.**